


X-Biotech™ Pre-stained Protein Ladder

1.Product Basic Information

Item	Specification
Cat. No.	X10116
Product Size	250 μL × 2
MW Range	8—180 kDa
Band Number	10
Storage	4°C for 2 months -20°C for 2 years
Storage Buffer	20 mM Tris-PO4 (pH 7.5), 2 mM EDTA.2 Na, 4 M Urea, 2 mM DTT,1.5% (W/V) SDS, 15% (V/V) Glycerol, 0.1% (V/V) Proclin 300

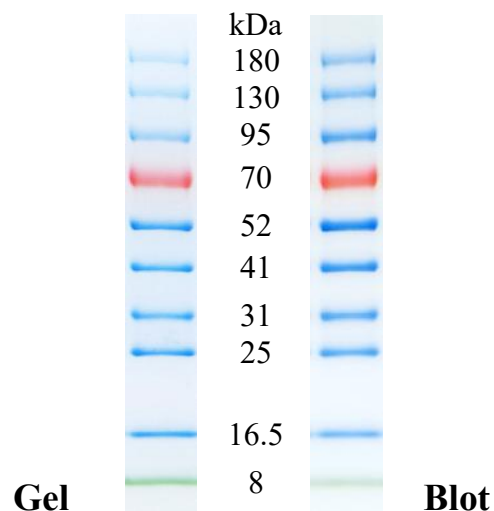

4-20% Tris-Glycine SDS-PAGE

Migration Patterns and approximate MWs (kDa) of X-Biotech™ Protein Ladder

Gel type	Tris-Glycine				Bis-Tris				Tris-Acetate		
	8%	10%	12.5%	15%	4-20%	4-12%	8-16%	4-20%	7%	3-8%	
Gel con- centration											
Running buffer	Tris-Glycine				MOPS				MES		Tris-Acetate
Apparent Molecular Weights, kDa											
10		180	180	180	180	180	180	180	180		
		130	130	130	130	130	130	130	130		
20	180	180	95	79	130	130	95	95	95		
	130	72	75	52	95	95	65	65	65		
30	130	52	41	41	70	65	52	52	52	180	
	95	41	31	31	52	52	41	41	41	130	
40	95	41	31	31	41	41	31	31	31	95	
	70	31	25	25	31	31	25	25	25	65	
50	70	31	25	25	25	25	16.5	16.5	16.5	180	
	52	25	16.5	16.5	16.5	16.5	8	8	8	130	
60	52	25	16.5	16.5	16.5	16.5	8	8	8	95	
	41	16.5	8	8	8	8	8	8	8	65	
70	41	16.5	8	8	8	8	8	8	8	52	
	25	8	8	8	8	8	8	8	8	41	
80	25	8	8	8	8	8	8	8	8	31	
	16.5	8	8	8	8	8	8	8	8	25	
90	16.5	8	8	8	8	8	8	8	8	16.5	
	8	8	8	8	8	8	8	8	8	8	
100	8	8	8	8	8	8	8	8	8	8	
	8	8	8	8	8	8	8	8	8	8	

2. Product Basic Information

The X-biotech™ Pre-stained Protein Ladder is a ready to use pre-stained mixture of 10 recombinant proteins ranging from 8 kDa to 180 kDa. It features the following key characteristics:

• Pre-stained Design

Three different chromophores are bound to the proteins, producing a brightly colored ladder. When separated by SDS-PAGE, most proteins exhibit a blue color, except for two reference bands: **70 kDa (orange-red)** and **8 kDa (green)**.

• Ready-to-Use Convenience

No additional steps (e.g., heating, dilution, or adding reducing agents) are required before use.

• Calibration Basis

The indicated apparent molecular weights are calibrated against unstained protein ladders from standard manufacturers, including *X-Biotech™ Protein Ladder (Cat. No. X10139)*, *Bio-Rad Laboratories (Cat. No. 1610363)* and *Thermo Fisher Scientific (Cat. No. 26614)*.

• Core Applications

1. Monitoring protein separation processes during SDS-PAGE electrophoresis.
2. Verifying Western blot transfer efficiency on membranes (PVDF or nitrocellulose membranes).
3. Approximating the molecular weight of target proteins on gels or blots.

• Shipping Condition

Transport with ice bag (4-8°C).

Note: Upon receipt of the product, please store it at -20°C promptly.

3. Recommended Experimental Procedures

3.1 Thawing and Preparation

Thaw the ladder at room temperature until completely dissolved (to redissolve any precipitated solids).

Gently but thoroughly mix the solution to ensure homogeneity.

Critical Warning: Do not heat the product above room temperature.

3.2 Recommended loading volumes

When performing polyacrylamide gel electrophoresis (SDS-PAGE), load an appropriate volumes of the ladder on an SDS-polyacrylamide gel:

- Mini gel electrophoresis: 5 µL per well
- Large gel electrophoresis: 10 µL per well

Use the same volumes for Western blotting.

The loading volumes listed above are recommended for gels with a thickness of 0.75-1.0 mm.

The loading volume should be doubled for 1.5 mm thick gels.

4. Key Notes and Precautions

4.1 Electrophoresis-Related Notes

1. In gels with a concentration <10%, low-molecular weight proteins in the ladder may migrate with the dye front.
2. Western Blot for Large Proteins (>100 kDa):
 - Extend transfer time or increase transfer voltage to ensure complete transfer.
 - SDS in Transfer Buffer: It is recommended to avoid adding SDS. If necessary for the experiment, the concentration must not exceed 0.02-0.04%.
3. Pre-stained proteins may show different apparent molecular weights in different SDS-PAGE buffer systems. For accurate approximate molecular weight determination:
 - Calibrate with unstained proteins in the same buffer system in advance.
 - Refer to the attached "Mobility Table" for migration patterns under different electrophoresis conditions.

4.2 Usage Restriction

For Research Use Only: This product is strictly for laboratory research purposes and not approved for diagnostic procedures (e.g., clinical sample testing, disease diagnosis).

5. Migration Patterns and Apparent Molecular Weights (kDa)

The apparent molecular weights under different gel types, gel concentrations, and running buffer conditions can be referenced in the "Migration Patterns and approximate MWs (kDa) of X-Biotech™ Protein Ladder".